


# MEASURING PICOSECOND FLUORESCENCE LIFETIMES

Using a Chromacity 520 in line with an Edinburgh Instruments FLS1000 Equipped with a Hybrid Photodetector

## Introduction

The measurement of fluorescence lifetimes provides a wealth of information on the sample under study, from charge-carrier lifetimes in semiconductors to the local environment of biomolecules in living cells. In many of these studies, short fluorescence lifetimes on a picosecond timescale are required to be measured.


In this application note, the measurement of sub 20 picosecond lifetimes is achieved by combining a Chromacity 520<sup>2</sup>, ultrashort pulse laser source with an Edinburgh Instruments FLS1000 Photoluminescence Spectrometer equipped with a hybrid photodetector.

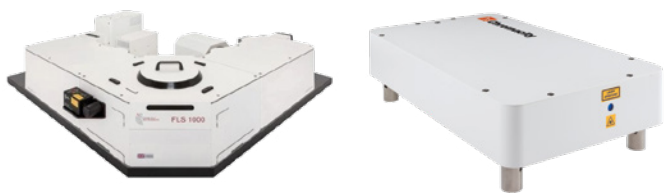

**FIGURE 1.** The Chromacity 520 was coupled into the FLS1000 Photoluminescence Spectrometer to narrow the IRF and drive performance.

## Materials

The hemicyanine dye trans-4-[4-(dimethylamino)-styryl]-1-methyl-pyridiniumiodide (4-DASPI) was chosen to demonstrate the measurement of short lifetimes using the FLS1000. DASPI dyes are commonly used as short lifetime fluorescence probes and as a laser gain media. 4-DASPI was purchased from Sigma Aldrich® and dissolved in either anhydrous ethanol or distilled water.

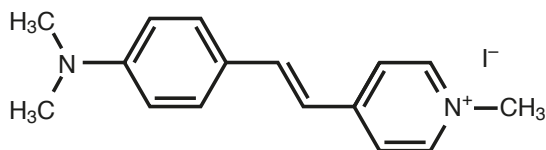

**FIGURE 2.** Molecular structure of 4-DASPI

## Experimental Configuration

Fluorescent lifetimes in the FLS1000 are measured using the time-correlated single photon counting (TCSPC) technique. In an ideal TCSPC system, the temporal width of the excitation pulse and the temporal response of the detection system would both be infinitely sharp (delta functions). However, this is never the case since lasers have finite pulse widths and detectors have transit response times. The deviation from the ideal is characterised by the instrument response function (IRF) of the TCSPC system. The IRF can be thought of as temporal response that the TCSPC system records for an infinitely short fluorescence lifetime.

The measured fluorescence decay is convoluted with the IRF, and the Full Width Half Maximum also known as the temporal width of IRF ( $FWHM_{IRF}$ ) determines the minimum lifetime that can be successfully measured.

A useful rule of thumb is that the minimum fluorescence lifetime that can be determined through reconvolution analysis is around 1/10th of the  $FWHM_{IRF}$ . It is therefore important to minimise the  $FWHM_{IRF}$  when measuring short fluorescence lifetimes.

The FLS1000 is a fully modular spectrometer and can be configured to suit different research requirements. By changing the configuration of the FLS1000 or the light source, the width of the IRF can be adjusted to meet the minimum measurable fluorescence lifetime required in different research applications.

**In this application note, the measurement of sub 20 picosecond lifetimes is achieved by combining a chromacity 520, ultrashort pulse laser source with an Edinburgh Instruments fls1000 photoluminescence spectrometer equipped with a hybrid photodetector.**

# MEASURING PICOSECOND FLUORESCENCE LIFETIMES

Using a Chromacity 520 in line with an Edinburgh Instruments FLS1000 Equipped with a Hybrid Photodetector

$$FWHM_{IRF} = \sqrt{FWHM_{detector}^2 + FWHM_{laser}^2 + FWHM_{dispersion}^3 + FWHM_{electronics}^2}$$

**FIGURE 3.** There are four main components which combine to give the total width of the IRF.


## FWHM Electronics

$FWHM_{electronics}$  is the broadening of the IRF due to the electronic jitter of the TCSPC timing electronics. This contribution to the IRF is small as every FLS1000 uses the advanced Edinburgh Instruments TCC2 counting electronics module which has a jitter as low as 20 ps.

## FWHM Dispersion

$FWHM_{dispersion}$  is the temporal dispersion of the light after passing through the emission monochromator. The extent of the dispersion depends on the groove density of the monochromator grating and the width of the monochromator slits. The higher the groove density the larger the spectral and temporal dispersion. High groove density gratings are therefore optimal for high resolution spectral measurements but broaden the IRF.

Each monochromator in the FLS1000 can house up to three gratings simultaneously enabling an optional low groove density grating to be installed for measuring short lifetimes. The user can quickly switch between the spectral (typically 1200 gr/mm) and lifetime (300 gr/mm) gratings in the Fluoracle® operating software.

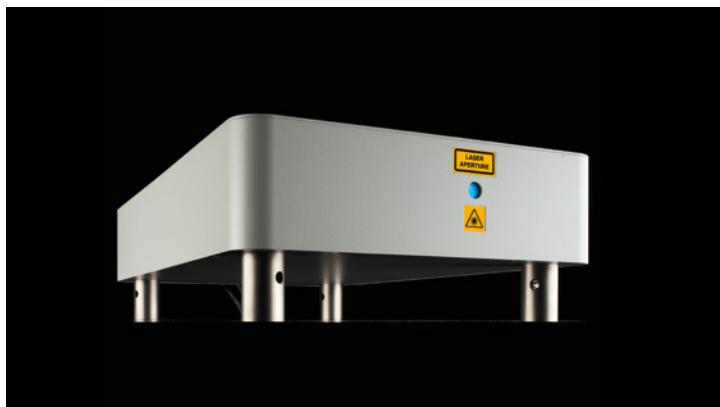

**FIGURE 4.** The Chromacity 520 is a compact, air cooled laser source making it easy to set up and use

For the measurement of 4-DASPI in the following section, the 300 gr/mm lifetime grating was used. To lower the temporal dispersion even further a subtractive double monochromator could be used, as found in the Lifespec II Spectrometer, which has zero temporal dispersion.

## FWHM Laser

$FWHM_{laser}$  is the pulse width of the laser excitation source. Picosecond pulsed diode lasers are the most common light source for TCSPC measurements but are not ideal for all scenarios. For that reason the FLS1000 has laser ports to allow the user to couple the Chromacity 520 into the spectrometer. By making use of this ultrashort laser source it has been possible to narrow the IRF. In this instance the Chromacity 520 generated pulse durations of 150 fs for the measurement of 4-DASPI. To trigger the TCSPC electronics, a fraction of the laser pulse was picked off and directed into the OT900 Optical Trigger Module.

## FWHM Detector

$FWHM_{detector}$  is broadening due to the transit time spread (TTS) of the detector and has the greatest influence on the overall width of the IRF.

The TTS is due to the photon-to-photon variation in the delay between the absorption of a photon at the photocathode of the detector and the electrical output pulse. The standard detector of the FLS1000 is the PMT-900 which is a conventional side window photomultiplier tube (PMT) which has a TTS of ~600 ps and limits the minimum lifetime to > ~60 ps. To measure shorter lifetimes the FLS1000 can be equipped with high-speed PMTs (HS-PMT) which have TTS of ~180 ps. For many years the only option for measuring shorter fluorescence lifetimes was using a microchannel plate PMT (MCP-PMT) which have excellent TTS of <25 ps but are expensive and fragile. Recently a new type of fast detector has been developed, the hybrid photodetector (HPD), which combines components of a traditional PMT with an avalanche photodiode to make a robust detector with a TTS as low as an MCP-PMT. For the measurement of 4-DASPI, the FLS1000 was equipped with an HPD detector. The combination of HPD detector, low groove density grating and 150 fs pulses from the Chromacity 520 laser resulted in a  $FWHM_{IRF}$  of only 34 ps.


# MEASURING PICOSECOND FLUORESCENCE LIFETIMES

Using a Chromacity 520 in line with an Edinburgh Instruments FLS1000 Equipped with a Hybrid Photodetector

## Results

The fluorescence decay of 4-DASPI in ethanol is shown in Figure 5a. The decay was measured at the emission maximum (600 nm) and the IRF was measured at the laser wavelength (520 nm) using a scattering dispersion of colloidal silica (LUDOX). The decay was fit with a single exponential using reconvolution fitting in the Edinburgh Instruments FAST advanced lifetime analysis software. A single exponential decay models the fluorescence well and the lifetime of 57 ps is in good agreement to previous reports.<sup>1</sup>

It is known that when 4-DASPI is dissolved in water its fluorescence lifetime decreases fivefold,<sup>1</sup> and a solution of 4-DASPI in water was therefore also measured to provide a more challenging sample to test the short lifetime capabilities of the FLS1000. The fluorescence decay and fit of 4-DASPI in water are shown in Figure 5b revealing a lifetime of 11 ps which is in perfect agreement to the 11 ps lifetime previously reported in the literature.<sup>3</sup>

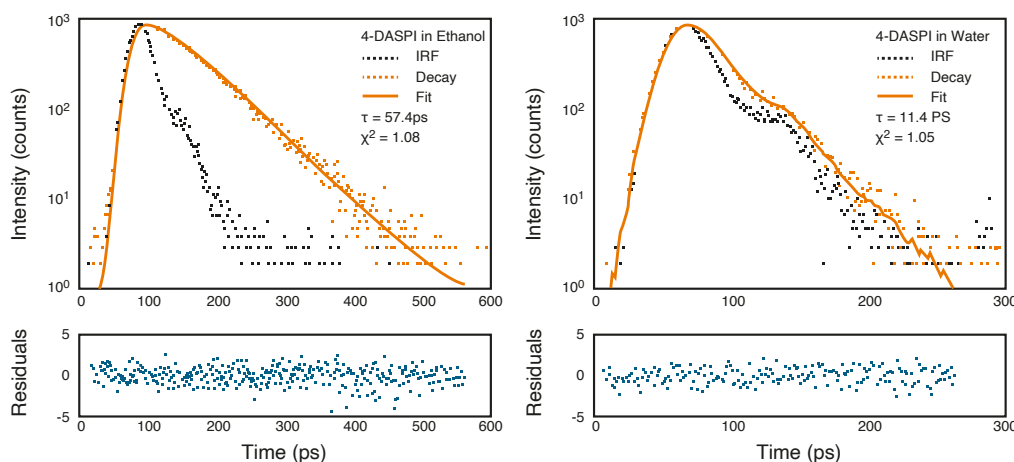

**FIGURE 5.** Fluorescence decays of 4-DASPI in (a) ethanol and (b) water measured using TCSPC.

## Conclusion

The capability of the FLS1000 Photoluminescence Spectrometer, in line with the Chromacity 520 ultrashort pulse laser, to measure short picosecond lifetimes was demonstrated by measuring the fluorescence lifetime of 4-DASPI in ethanol (57 ps) and water (11 ps). Through customisation of the light sources, gratings and detectors of the modular FLS1000, the minimum measurable lifetime can be adjusted to meet the demands of a wide variety of different research areas.

*Content courtesy of Edinburgh Instruments*

## References.

1. J. Kim, M. Lee, J.-H. Yang, & J.-H. Choy, Photophysical Properties of Hemicyanine Dyes Intercalated in Na-Fluorine Mica, *J. Phys. Chem. A* 104 1388-1392 (2000)
2. Chromacity 520 Data Sheet, [www.chromacitylasers.com](http://www.chromacitylasers.com)
3. J. Kim & M. Lee, Excited-State Photophysics and Dynamics of a Hemicyanine Dye in AOT Reverse Micelles, *J. Phys. Chem. A* 103 3378-3382 (1999)

Learn how our ultrafast lasers can enable you to discover more.  
For more information, email: [sales@chromacitylasers.com](mailto:sales@chromacitylasers.com)

